


pipetman®

DIAMOND TIPS

Refill for TOWERPACK boxes

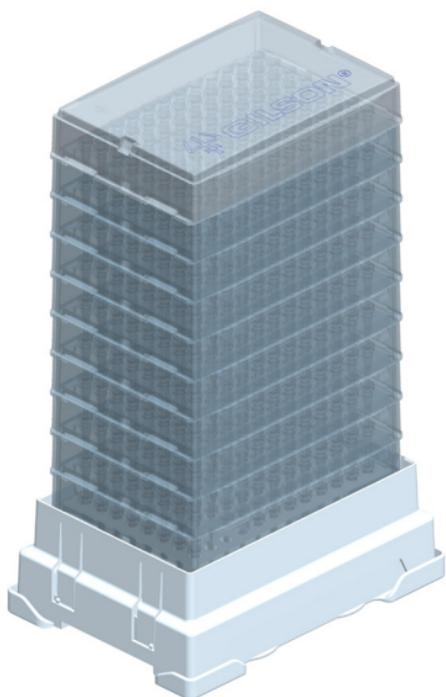

- Uses 60% less space and 70% less plastic
- One reload box for the entire TOWERPACK range
- Easy-to-use, one-handed system
- Protective base to prevent contamination from environmental pollution and dust
- Available in standard and sterilized versions

Loading instructions on the reverse side.



Made in France

Download certificates at www.gilson.com


Gilson S.A.S.

19, avenue des Entrepreneurs BP 145 • F-95400
Villiers-le-Bel, France
T +33 (0) 1 34 29 50 00 • F +33 (0) 1 34 29 50 20

 **GILSON®**

Loading a TOWERPACK Box

1. Remove the TOWERPACK dispenser from the protective support.
2. Place the TOWERPACK on the opened reload box.
3. Press down firmly until the rack snaps into place.
4. When a rack is empty, remove it from the reload box, and load a new rack.

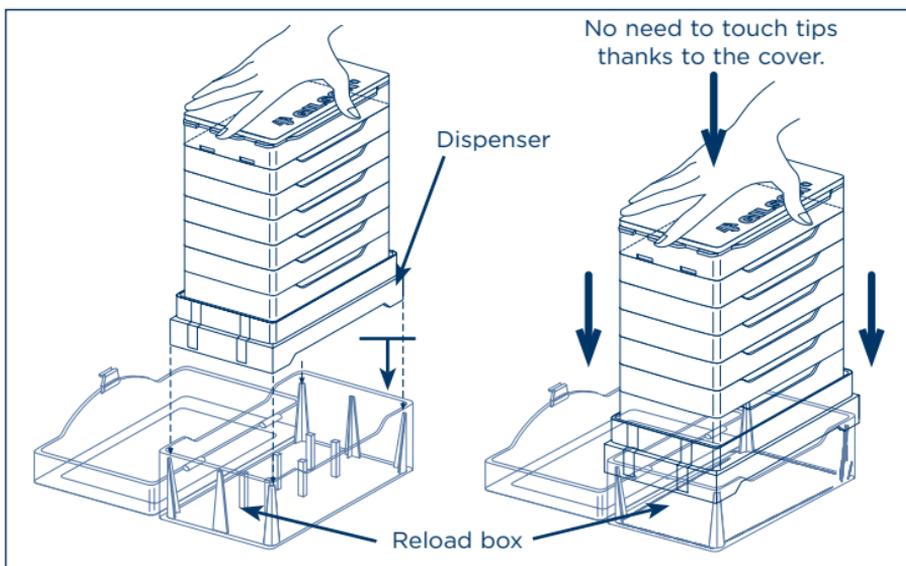

Autoclaving

The entire TOWERPACK can be autoclaved or just the reload box filled with one rack of tips.

Autoclaving conditions: moist heat, 121°C, 0.1 MPa for 20 minutes.

NOTE

When autoclaving PIPETMAN® DIAMOND Tips D10, keep the box upright to prevent these tips from falling out of the rack.

Loading a Multichannel PIPETMAN®

The ROCKY RACK technique makes it extremely easy to fit PIPETMAN DIAMOND Tips on a multichannel PIPETMAN.

Exert a light vertical force to obtain a perfect seal on all eight or twelve channels.



ROCKY RACK technique